

GET THE MOST OUT OF YOUR MICROSCOPE

Upgrade your ZAP with new features and apps

Here's why you should upgrade – and how.



ACTIVATE NEW FEATURES INSTEAD OF BUYING A NEW MICROSCOPE

Why you should upgrade

Perhaps your inspection processes have changed and you need help maintaining your product quality and efficiency. Or maybe you've realized there's more to digital microscopy than you first thought. In either case, here's why you should upgrade your ZAP to a MOVE instead of purchasing a new microscope.

If you want to maintain the flexibility of the ZAP but need the advanced features, apps, presets and documentation options of the TAGARNO TREND and PRESTIGE, you can purchase a MOVE upgrade kit.

The MOVE Upgrade kit will transform the ZAP microscope into an advanced and customizable microscope – at a fraction of the price compared to purchasing a new microscope.

With the MOVE upgrade kit and the included Ruler app, Focus stacking app and Watermark app, you'll be able to:

- Make on-screen measurements
- Create ultra-sharp photos with no blurred or out of focus areas
- Save photos with your logo on

You will also gain access to unprecedented processing power that is guaranteed to take your inspections to the next level.

There's two upgrade processes depending on when you bought your ZAP.

Read more about that on the next pages.

SCENARIO 1

You're ready to go in 3 simple steps

If your microscope meets one of the following criteria, all you need to upgrade from ZAP to MOVE is a MOVE upgrade kit:

- Has serial number 33504 or higher*
- Has Firmware version 5.4

STEP 1

Buy a MOVE Upgrade kit**

STEP 2

Connect it to your ZAP

STEP 3

Use your TAGARNO
MOVE


** The serial number is available in the top left corner on the monitor when the microscope is powered on as well as on the back of the camera head*

*** Ask your distributor for a quote on the MOVE Upgrade kit*

SCENARIO 2

A bit of help is needed before you can use your MOVE

If your microscope does not run on Firmware version 5.4 or has a serial number below 33504*, a spare part in the microscope needs to be changed before you can get the benefits of a MOVE.

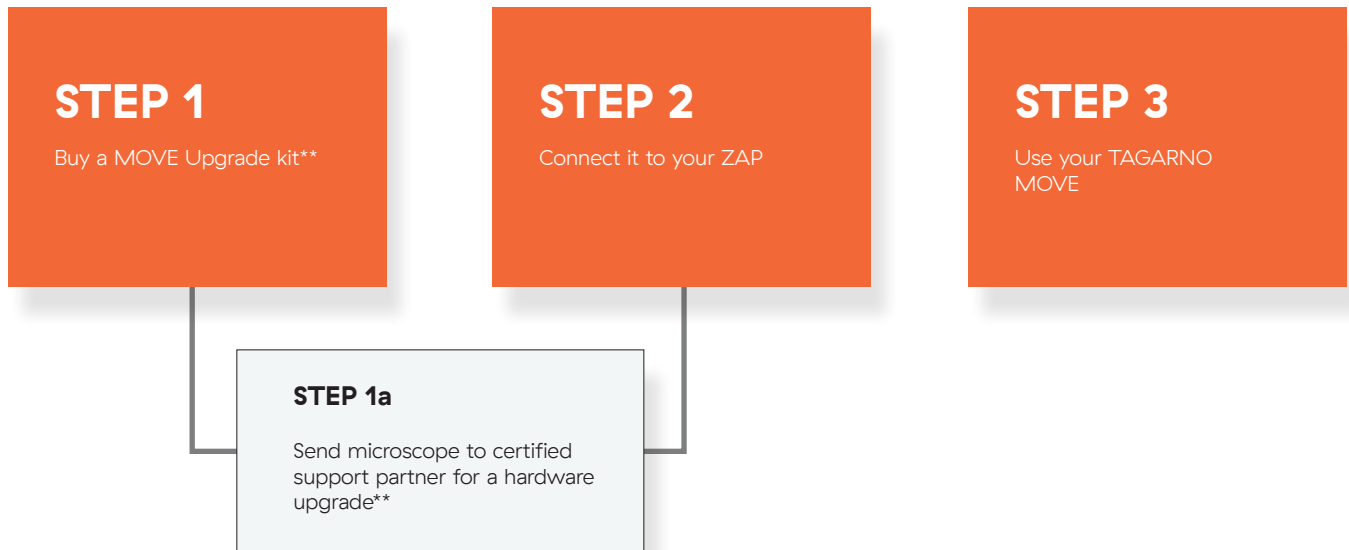

* The serial number is available in the top left corner on the monitor when the microscope is powered on as well as on the back of the camera head

** Ask your distributor for a quote on the MOVE Upgrade kit and hardware replacement

WANT TO KNOW MORE?

Contact us for more info

TAGARNO MOVE is the fastest and cheapest way to optimize your inspection processes, regardless of the upgrade process you need.

Visit tagarno.com or contact your local TAGARNO distributor if you have any questions on the TAGARNO MOVE. We're here to help!

TAGARNO